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TEL: 0748-20-3535 FAX: 0748-25-7171

Mail: biosales@atect.co.jp

atect Stackable Medium for aerobic count "aS-Medium AC" instruction manual

Product Information

aS-Medium AC consists of a plastic substrate and a cover film. As an alternative to the conventional method using an agar medium, it is an easy to use microbiological test medium that measures the aerobic count from a sample solution. Plates can be stacked stably due to the indentations at the four corners. In addition, since this product is in the form of a thin sheet, it contributes to space saving during storage and reduction of environmental load during disposal.

aS-Medium AC is intended to be used by trained laboratory specialists or by general users in the presence of those specialists in a microbiology laboratory setting.

Materials Required, not Provided

1. Incubator (35±1°C).
2. Stomacher.
3. Stomacher bag with filter (250 or 500 mL volume).
4. Butterfield's phosphate-buffered dilution water
5. Pipettor, pipette, and tips

Sample Preparation

1. Weigh 50 g of the sample and add it to the filtered Stomacher bag.
2. Add 450mL of the appropriate buffer solution (such as Butterfield's phosphate-buffered diluent) to the sample to filtered Stomacher bag.
3. Use a Stomacher to homogenize the sample.
※Alternatively, for steps 1 to 3 above, 2 x 25 g portions can be weighed into separate stomacher bags. Add 225 mL Butterfield's phosphate-buffered diluent to each portion and stomach for 2 min. After homogenization, combine the two portions into a single container and briefly mix by hand.
4. Dilute the sample solution as needed.

Inoculation

1. Lift the cover film and pipette 1 mL of the sample solution onto the medium. Return the cover film to cover the medium surface.
2. Let stand for 3 minutes.
3. Stack the plates with the cover film on top.
It can be stacked up to 25 plates.
4. Incubate the plate at 35±1°C for 48±2h.

Interpretation

The surface of the cover film facing the medium contains the color-developing agent tetrazolium salt (2,3,5-Triphenyl Tetrazolium Chloride). Count the colonies that have turned red.

Precautions for counting colonies

The area of the medium is about 20 cm². The range of the number of bacteria that can be counted is 30 to 300 cfu. If the number of colonies is large, it can be estimated by counting 1 square cm (1 cm x 1 cm) in the grid and multiplying by 20.

Storage and Shelf life

1. Store the product in a refrigerator at 4 - 10°C.
2. Shelf life is 3 months (stored at 4 - 10°C, unopened).
3. The expiration date is written on the lower left of the product in addition to the label on the aluminum bag.
4. After opening, return it to the aluminum bag and store it in the refrigerator (4 - 10°C) in a sealed state, and use it within 1 month.

Precautions for use

1. Check the instruction manual before using this product.
2. This product is mainly used for microbiological testing of food and the environment samples. Do not use it for purposes other than the intended purpose.
3. The product should only be handled by experienced individuals.
4. This product has not been verified for all foods, inspection procedures, and strains.
5. Please note that the user is responsible for the interpretation and operation of the inspection results, and the manufacturer is not responsible for it.
6. Products that have passed the expiration date are not covered by the warranty.
7. If the product is damaged, foreign matter, discolored, etc., refrain from using it and contact the manufacturer / distributor.

Safety Precautions

1. Assay users should observe standard microbiological practices and safety precautions when performing this assay.
2. Should product or reagents get in your eyes or mouth, rinse thoroughly with water and consult a doctor.
3. Used medium must be disposed after autoclaving. It is the responsibility of the end user to handle waste and effluents processed according to their nature and degree of hazard and to treat and dispose of them in accordance with applicable local, state, and federal regulations.

Manufacturer information

Atect Corp.

Address: 3275-1, Kamihaneda-cho, Higashiomi city, Shiga, 527-0082, Japan

TEL: 0748-20-3535 FAX: 0748-25-7171

Mail: biosales@atect.co.jp