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TEL: 0748-20-3535 FAX: 0748-25-7171

Mail: biosales@atect.co.jp

## atect Stackable Medium for aerobic count "aS-Medium AC" instruction manual

### Product Information

aS-Medium AC consists of a plastic substrate and a cover film with a resin frame. As an alternative to the conventional method using an agar medium, it is an easy to use microbiological test medium that measures the aerobic count from a sample solution. Plates can be stacked stably due to the indentations at the four corners. By adding an adhesive resin frame to the cover film, it improves adhesion to the substrate, prevents the cover film from unintentionally turning over, and suppresses leakage of the test liquid. In addition, since this product is in the form of a thin sheet, it contributes to space saving during storage and reduction of environmental load during disposal.

aS-Medium AC is intended to be used by trained laboratory specialists or by general users in the presence of those specialists in a microbiology laboratory setting.

### PTM Certification

aS-Medium AC has been evaluated against a reference method (MLG3.02) in the AOAC Research Institute PTM program and is certificated as AOAC PTM 112202 for enumeration of mesophilic aerobic bacteria in meat products (ready-to-eat ham, ready-to-eat bacon, fresh raw pork, fresh raw beef trim, and fresh raw ground beef). For evaluation in this test, 2 sheets are used for each test area, and the number of bacteria is calculated based on the average value and compared with the reference method.

### Materials Required, not Provided

1. Incubator ( $35\pm1^{\circ}\text{C}$ ).
2. Stomacher.
3. Stomacher bag with filter (250 mL or 500 mL volume).
4. Butterfield's phosphate-buffered dilution water.
5. Pipettor, pipette, and tips.

### Sample Preparation

1. Weigh 50 g of the sample and add it to the filtered Stomacher bag.
2. Add 450mL of the appropriate buffer solution

(such as Butterfield's phosphate-buffered diluent) to the sample to filtered Stomacher bag.

3. Use a Stomacher to homogenize the sample.  
※Alternatively, for steps 1 to 3 above, 2 x 25 g portions can be weighed into separate stomacher bags. Add 225 mL Butterfield's phosphate-buffered diluent to each portion and stomach for 2 min. After homogenization, combine the two portions into a single container and briefly mix by hand.
4. Dilute the sample solution as needed.

### Inoculation

1. Pinch the cover film between your thumb and index finger and place your remaining fingers from the center of the film over the logo in the upper left corner.
2. Use your thumb and index finger to turn over the bottom edge of the resin frame of the cover film (the adhesive strength at the bottom edge of the resin frame is strong, so pinch the film firmly and apply force to turn it over).
3. Turn over the cover film while holding down the product. (Do not remove the cover film completely. Leave the top edge glued.)
4. Drop 1 mL of the sample solution onto the medium by pipette. Return the cover film to cover the medium surface.
5. Let stand for 3 minutes.
6. Stack the plates with the cover film on top.  
It can be stacked up to 25 plates.
7. Incubate the plate at  $35\pm1^{\circ}\text{C}$  for  $48\pm2\text{h}$ .

### Interpretation

The surface of the cover film facing the medium contains the color-developing agent tetrazolium salt (2,3,5-Triphenyl Tetrazolium Chloride). Count the colonies that have turned red.

Conduct the test using 2 sheets per test area based on the test method used for PTM certification. Calculate the average number of bacteria on each plate and calculate the number of bacteria in the sample from that number.

### Precautions for counting colonies

The area of the medium is about 20 cm<sup>2</sup>. The range of the number of bacteria that can be counted is 30 to 300 cfu. If the number of colonies is large, it can be estimated by counting 1 square cm (1 cm x 1 cm) in the grid and multiplying by 20.

### Storage and Shelf life

1. Store the product in a refrigerator at 4 – 10°C.
2. Shelf life is 13 months (stored at 4 – 10°C, unopened).
3. The expiration date is written on the lower left of the product in addition to the label on the aluminum bag.
4. After opening, return it to the aluminum bag and store it in the refrigerator (4 - 10°C) in a sealed state, and use it within 1 month.

### Precautions for use

1. Check the instruction manual before using this product.
2. This product is mainly used for microbiological testing of food and environmental samples. Do not use it for purposes other than the intended purpose.
3. The product should only be handled by experienced individuals.
4. Do not place the product under an ultraviolet lamp after taking it out of the aluminum bag. The ingredients applied to the cover film may react and turn red. There is no problem using it under a fluorescent light.
5. This product has not been verified for all foods, inspection procedures, and strains.
6. Some bacterial strains, such as lactic acid bacteria and Micrococcus, may have low detection sensitivity or may not be detected. Additionally, for some strains, higher numbers of bacteria may be detected than with plate count agar.
7. Please note that the user is responsible for the interpretation and operation of the inspection results, and the manufacturer is not responsible for it.
8. Products that have passed the expiration date are not covered by the warranty.
9. If the product is damaged, foreign matter, discolored, etc., refrain from using it and contact the manufacturer / distributor.

### Safety Precautions

1. Assay users should observe standard microbiological practices and safety precautions when performing this assay.
2. Should product or reagents get in your eyes or mouth, rinse thoroughly with water and consult a doctor.
3. Used medium must be disposed after autoclaving. It is the responsibility of the end user to handle waste and effluents processed according to their nature and degree of hazard and to treat and dispose of them in accordance with applicable local, state, and federal regulations.

### Manufacturer information

Atect Corp.

Address: 3275-1, Kamihaneda-cho, Higashiomi city, Shiga,

527-0082, Japan

TEL: 0748-20-3535 FAX: 0748-25-7171

Mail: biosales@atect.co.jp